

Antibody Pairing Results for CPTC-PDLIM1

| “Capture” (bottom) Antibody | “Indicator” (top) Antibody | | |
|--------------------------------|----------------------------|----------|----------|
| | PDLIM1-1 | PDLIM1-2 | PDLIM1-3 |
| PDLIM1-1 | - | ++++ | ++++ |
| PDLIM1-2 | + | - | +++ |
| PDLIM1-3 | ++++ | + | - |

Method – BioRad ProteOn XPR:

- Surface coated with Rabbit anti-mouse IgG Fc specific antibody
- “Bottom” antibody flowed over the surface
- Antigen flowed over the surface
- Orthogonal flow of “Top” antibody over the surface

Results:

- No pairing observed
- + Weak Pairing
- ++ Moderate Pairing
- +++ Strong pairing